


# Product Data Sheet

Hand refractometer  
Reference 95000-015

## Specifications


- 0-90 °Brix
- Div 0.5°Brix
- Classic metal - blue line
- In box
- With cloth
- Pipette
- Pocket case
- Screwdriver
- Instructions for use

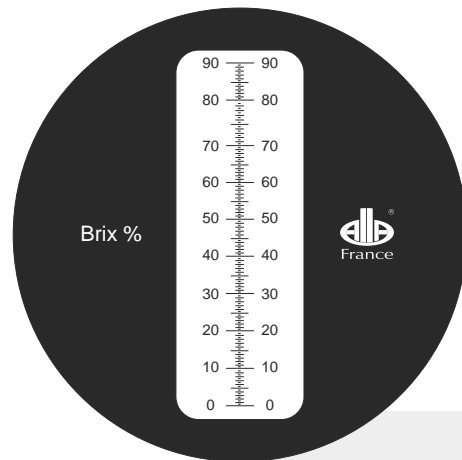

## Instructions

### Components

- 1- Prism
- 2- Cover plate
- 3- Setting screw
- 4- Mirror tube
- 5- Eyepiece (adjusting ring of dioptre)

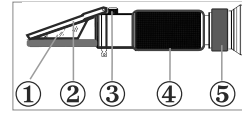

### Calibration

Aim the front end of the refractometer in the direction of bright light and set the adjusting ring of dioptre until the scale can be seen clearly. Open the cover plate, put one or two drops of standard liquid on the prism. Close the cover plate and press it lightly. Then adjust the setting screw to make the light and dark limit coincide with the 60°Brix point. If the instrument is ATC (with Automatic Temperature Compensation system), the room temperature must be 20°C whenever the instrument is recalibrated. Once calibrated, shifts in ambient temperature within the acceptable range (between 10°C/30°C) should not effect accuracy. The calibration of our refractometers can be made with our master liquids which are available for purchase.

### Utilisation

Open the cover plate, clean the surface of the prism with a soft cotton cloth, put one or two drops of the solution to be measured onto the prism. Close the cover plate and press it lightly, then read the corresponding scale of light and dark limit indicating the value of the measured solution. In case of the temperature has a lot of variation, you should control the reference point every 30 minutes. After the measurement, clean the remaining liquid on the surface of the prism and the cover plate with a damp gauze. After drying, store the instrument carefully.

### Automatic Temperature Compensation system

The reference temperature is 20°C (68°F). Refractometers with "ATC" have an automatic temperature compensation.

### Caution

Do not wash the instrument with water in order to avoid water entering inside. Handle gently and do not touch the optical surfaces. The refractometer must be stored in a dry, clean and non corrosive environment. Avoid strong shocks.

Alla France

Z.I. du Bompas - 13, rue du Bompas - Chemillé - 49120 Chemillé-en-Anjou - FRANCE  
Tél. + 33 (0)2 41 30 55 08 - Fax +33 (0)2 41 30 34 67

[www.allafrance.com](http://www.allafrance.com)

  
France